

4.4 Adjusting the Rocking Speed

Start the Rocker II by activating the power switch. To increase rocking speed adjust the speed control clockwise.

5. Accessories

5.1 Staking Tray 12" x 12" with flat mat

Boekel Part Number 260355

5.2 Dimpled Mat 12" x 12" (vinyl)

Boekel Part Number C2202208

5.3 Flat Mat 12" x 12" (rubber)

Boekel Part Number C2202162

6. Fault Diagnosis

Symptom

1. Unit does not operate

Possible Cause

- a. Unit not switched on
- b. Unit not plugged into power supply
- c. Power supply failure

Action Required

- a. Switch on
- b. Plug in, switch on
- c. Check that other electrical appliances on the same circuit are working

7. Technical Specifications

This equipment is intended for indoor use and will meet its performance figures within the ambient temperature range of +4°C to 40°C, with maximum relative humidity of 80% (non-condensing). Installation Category II (transient voltages). Pollution Degree 2 in accordance with IEC 664. Suitable for operation at altitudes of up to 6500 feet.

Specifications:

Temperature Range:	+4°C to 40°C
Supply Voltage Range:	115V +/- 10%, 50/60 Hz 230V +/- 10%, 50/60 Hz
Power Rating:	Model 260350: 22 W Model 260350-2: 22 W
Load Capacity (with Stacking Platform)	2.7 kilogram (6 lbs.) 2.3 kilogram (5 lbs.)